

ON Sequence

1. Main Switch Power ON (left, black table)
2. Polychrome V (shelf): wait for ~ 1-2 minutes clicking sounds
3. ICU (normal table, left, lower)
4. Power supply scanner (normal table, left, higher)
5. SCU (black table) wait for clicking sound
6. Andromeda (IF NEEDED)
7. Clara Camera
8. Lasers (table, big black box)
9. Computer
10. Log-in qNumber and qPassword
11. Start Omicron Control Software
12. Start LA Software (Offline Analysis will be started automatically)

OFF Sequence

1. Close Omicron Software
2. Close Offline Analysis Software
3. Move down the objective in LA software > enter z=0
4. Close LA Software
5. Shutdown Windows
6. SCU (black table)
7. Power supply scanner (normal table, left, higher)
8. ICU (normal table, left, lower)
9. Laser (table, big black box)
10. Andromeda
11. Clara Camera
12. Polychrome V (shelf): DON'T RESTART IT WITHIN five minutes
13. Main Switch Power ON (left, black table)